

TRACEABLE® MINI THERMOMETER INSTRUCTIONS

SPECIFICATIONS

Range: –58 to 302°F (–50 to 150°C)

Resolution 0.1° from –20.0 to 100°C

Accuracy: ±1°C from –20 to 200°C

OPERATION

1. Slide the power “ON/OFF” switch to the “ON” position.
2. Slide the “C/F” switch to select the desired reading.
3. Press “HOLD” button once to freeze the display.
4. Press the “HOLD” button again to resume measurement.
5. Slide the power “ON/OFF” switch to the “OFF” position to conserve battery life.
6. Keep plastic housing and plastic handle away from heat source. Do not place inside oven or expose to open flame.
7. Do not immerse probe handle or cable in liquid.

BATTERY REPLACEMENT

1. Use a coin to open the battery cover on back of unit.
2. Replace with catalog number 1039.

WARRANTY, SERVICE, OR RECALIBRATION

For warranty, service, or recalibration, contact:

CONTROL COMPANY

4455 Rex Road
Friendswood, Texas 77546 USA
Ph. 281-482-1714 Fax 281-482-9448
E-mail sales@control3.com
www.control3.com

Control Company is ISO 9001 Quality-Certified by DNV and ISO 17025
accredited as a Calibration Laboratory by A2LA.