

TRACEABLE® WATER-RESISTANT, FLASHING TIMER INSTRUCTIONS

SPECIFICATIONS:

Display: 1/2" LCD
Range: 99 Minutes, 59 seconds to zero
Resolution: 1 second
Accuracy: 0.01%
Alarm: 1 Minute audible alarm, bright red LED flashing visual alarm
Function: Countdown timing
Features: Water-resistant, Audible alarm, Flashing red LED when time is up, Infinite timeouts, Supplied with lanyard and suction cup.

CONTROLS/INDICATORS

Separate START and STOP buttons, and ten time-setting buttons, numbered "1" through "10".

OPERATION

1. Select desired countdown time with buttons "1" through "10".
2. Press the "START" button to begin countdown.
3. Press the "STOP" button for timeouts or to stop the countdown.
4. When zero is reached a bright red LED flashes and an audible alarm sounds for one minute.
5. Alarm can be turned off by pressing any of the buttons on the unit.
6. To clear the time entry, press the "STOP" button twice.

ALL OPERATIONAL DIFFICULTIES

If this unit does not function properly for any reason, replace the battery with a new high-quality battery (see "Battery Replacement" section). Low battery power can occasionally cause any number of "apparent" operational difficulties. Replacing the battery with a new fresh battery will solve most difficulties.

BATTERY REPLACEMENT

Erratic readings, a faint display or no display are all indicators that the batteries must be replaced. The two battery compartments located on the back of the unit can be opened with a screw driver (turn in the direction indicated). Replace exhausted batteries with two new 1.5 volt batteries. When replacing battery compartment covers, carefully ensure that each cover is properly in place in order to maintain the unit's water-resistance. Replacement Battery Cat. No. 1039.

WARRANTY, SERVICE, OR RECALIBRATION

For warranty, service, or recalibration, contact:

CONTROL COMPANY

4455 Rex Road
Friendswood, Texas 77546 USA
Ph. 281-482-1714 Fax 281-482-9448
E-mail sales@control3.com
www.control3.com

Control Company is ISO 9001 Quality-Certified by DNV
and ISO 17025 accredited as a Calibration Laboratory by A2LA.

Cat. No. 5036

Traceable® Water-Resistant Flashing Timer Instructions

92-5036-00

Rev. 1 062504

Traceable® is a registered trademark of Control Company

© 2004  Control Company