


## INSTRUCTIONS FOR ASSEMBLY

### MODEL C-10 SERIES CHROMATO-VUE® CABINET

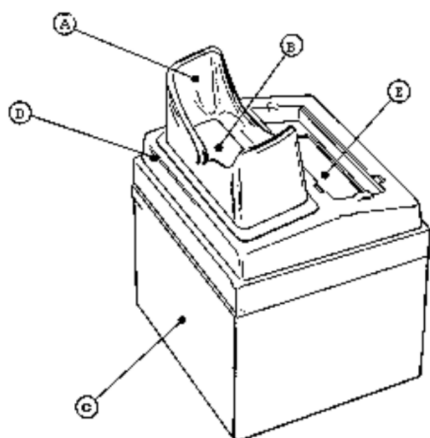

1. Lift flexible eyepiece (A) and cover assembly (D) out through top of cabinet.
2. Slip lamp well (E) portion of cover assembly over rim at back of cabinet until flush with top. Apply enough pressure toward front portion to snap over rim. Ensure that cover assembly is flush with top on all sides.
3. Slip flexible eyepiece over rim above contrast control filter (B) with notched portion facing front of cabinet.
4. Place the Mineralight® lamp into the molded lamp well with power cord placed to the left. The Chromato-Vue cabinet is now ready to use.
5. Lift light curtain (C) when placing or positioning materials; leave in closed position when in use to block external light from entering the cabinet.



For additional information on UVP's products, go to [www.UVP.com](http://www.UVP.com)

Corporate Headquarters: UVP, LLC  
2066 W. 11th Street Upland, CA 91786 / Tel: (800) 452-6788  
(909) 946-3197 / Fax: (909) 946-3597

European Operations: Ultra-Violet Products Ltd  
Unit 1, Trinity Hall Farm Estate, Nuffield Road,  
Cambridge CB4 1TG UK / Tel: +44(0)1223 420022  
Fax: +44(0)1223 420156



Replacement Parts:  
Eyepiece: 23-0036-01  
Contrast Control Filter: 38-0013-01

4. Place the portable lamp in the lamp support bracket. This bracket firmly holds the lamp in place. The Chromato-Vue cabinet is now ready to use. Lift light curtain (C) when placing or positioning materials; leave the curtain in the closed position when in use to block out any outside light.

To assemble the C-10P, follow the C-10 Cabinet Series instructions numbers 1 through 3 on the previous page. Instruction number 4 for the C-10P is as follows:

The difference between the C-10, C-10E and C-10P Cabinets is the addition of a lamp support bracket on the C-10P. The purpose of the bracket is to hold the portable lamp securely in place.

## C-10P CHROMATO-VUE CABINET

For additional information on UVP's products, go to [www.UVP.com](http://www.UVP.com)

Corporate Headquarters: UVP, LLC  
2066 W. 11th Street Upland, CA 91786 / Tel: (800) 452-6788  
(909) 946-3197 / Fax: (909) 946-3597

European Operations: Ultra-Violet Products Ltd  
Unit 1, Trinity Hall Farm Estate, Nuffield Road,  
Cambridge CB4 1TG UK / Tel: +44(0)1223 420022  
Fax: +44(0)1223 420156



Replacement Parts:  
Eyepiece: 23-0036-01  
Contrast Control Filter: 38-0013-01

4. Place the portable lamp in the lamp support bracket. This bracket firmly holds the lamp in place. The Chromato-Vue cabinet is now ready to use. Lift light curtain (C) when placing or positioning materials; leave the curtain in the closed position when in use to block out any outside light.

To assemble the C-10P, follow the C-10 Cabinet Series instructions numbers 1 through 3 on the previous page. Instruction number 4 for the C-10P is as follows:

The difference between the C-10, C-10E and C-10P Cabinets is the addition of a lamp support bracket on the C-10P. The purpose of the bracket is to hold the portable lamp securely in place.

## C-10P CHROMATO-VUE CABINET