

TRACEABLE® BIG-DIGIT 4-ALERT THERMOMETER °C/°F INSTRUCTIONS

OPERATING INSTRUCTIONS

Normal display shows inside and outside temp in 1/10 degree.



1. One point temp control


a. High Temp Alert "▲"

— Slide switch to temp "Alert" position, press "Alert Select" to select High Temp Alert "▲" mode.

- Press "Hi Set" to set desired High Temp Alert.
- Press "Alert On/Off" to turn on the High Temp Alert. ("Alert" flag appears on display.)
- Alarms and LCD flashes 5 sec every one min when the sensed temp is higher than the set temp.

b. Low Temp Alert "▼"

— Slide switch to temp "Alert" position, press "Alert Select" to select Low Temp Alert "▼" mode.

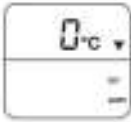

- Press "Lo set" to set desired Low Temp Alert.
- Press "Alert On/Off" to turn on the Low Temp Alert. (Alert flag appears on display.)
- Alarms and LCD flashes 5 sec every one min when the sensed temp is lower than the set temp.

FITTING THE OUTSIDE SENSOR: LOCATE OUTSIDE SENSOR AWAY FROM DIRECT SUNLIGHT AND HEAVY SATURATION.




2. Two point temp control

a. Inside Temp Alert "⌂"

— Slide switch to temp "Alert" position, press "Alert Select" to select Inside Temp Alert "⌂" mode.

- Press "Hi set" to set the desired upper temp alert value.
- Press "Lo set" to set the desired lower temp alert value.
- Press "Alert on/off" to turn, on the Inside Temp Alert. ("Alert" flag appears on display).
- Alarms and LED flashes 5 sec every one min when the sensed temp is within the set temp.

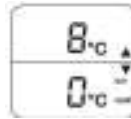

b. Outside Temp Alert "◆"

— Slide switch to temp "Alert" position, press "Alert select" to select Outside Temp Alert "◆" mode.

- Press "Hi set" to set the desired upper temp alert value.
- Press "Lo set" to set the desired lower temp alert value.
- Press "Alert on/off" to turn on the Outside Temp Alert. ("Alert" flag appears on display)
- Alarms and LED flashes 5 sec every one min when the sensed temp is outside the set temp.

3. Slide switch to "Temp" position after setting temp alert.

4. Alert off

Press "Alert on/off" in temp "Alert" position, "Alert" flag disappears on display.




Folding stand for counter top display

BATTERY REPLACEMENT:

- 1 Pc "AAA" or "UM4" battery
- Insert the battery as direction indicated.

Use fingernail to lift up battery cover and insert battery as to indicated direction.

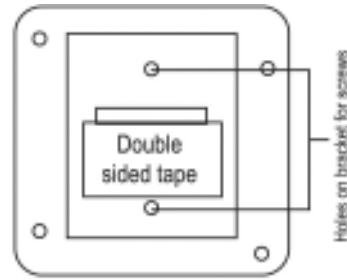

EASY MOUNTING OF BRACKET ON WALL BY:

- Double-sided tape or
- Screws

WARRANTY, SERVICE, OR RECALIBRATION

For warranty, service, or recalibration, contact:

CONTROL COMPANY

4455 Rex Road

Friendswood, Texas 77546 USA

Ph. 281-482-1714 Fax 281-482-9448

E-mail sales@control3.com

www.control3.com

Control Company is ISO 9001 Quality-Certified by DNV and ISO 17025 accredited as a Calibration Laboratory by A2LA.